




## Test Barcode (for Use with Scan Now)

### Instructions for Use

Please print this page\*. It includes an example Smart Label barcode with which you can test the Scan Now App and/or your USB barcode scanner.

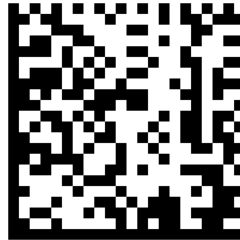


Example Barcode

**\*Note** – many mobile phone cameras and USB scanners can scan **directly from the screen**. To save paper you might want to try to scan from your screen before you print anything.

### Scan Now App Users

Download and launch the Scan Now app on your camera-enabled mobile device. Scan the test code by tapping the Scan Barcode button circled below:

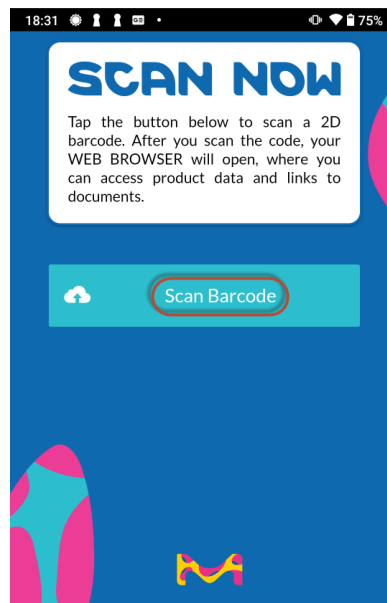

### USB Scanner Users

1. Go to the Scan Now page at <https://www.emdmillipore.com/ScanNow>
2. To scan your barcode, please use a standard barcode reader that supports GS1 two-dimensional (2D) barcodes. For best results, please set up your scanner to use an English keyboard layout.