




Product Information

SIGMA QUALITY CONTROL TEST PROCEDURE Enzymatic Assay of **SORBITOL DEHYDROGENASE** (EC 1.1.1.14)

PRINCIPLE:

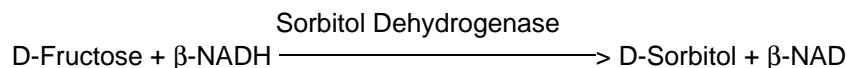

Abbreviations used:

β -NADH = β -Nicotinamide Adenine Dinucleotide, Reduced Form

β -NAD = β -Nicotinamide Adenine Dinucleotide, Oxidized Form

CONDITIONS: T = 25°C, pH 7.6, $A_{340\text{nm}}$, Light path = 1 cm

METHOD: Continuous Spectrophotometric Rate Determination

REAGENTS:

- A. 100 mM Triethanolamine Buffer, pH 7.6 at 25°C
(Prepare 100 ml in deionized water using Triethanolamine, Hydrochloride, Sigma Prod. No. T-1502. Adjust to pH 7.6 at 25°C with 5 M NaOH.)
- B. 1.1 M D-Fructose Solution (Fructose)
(Prepare 5 ml in deionized water using D(-)Fructose, Sigma Prod. No. F-0127.)
- C. 12.8 mM β -Nicotinamide Adenine Dinucleotide, Reduced Form, Solution (β -NADH)
(Dissolve the contents of one 10 mg vial of β -Nicotinamide Adenine Dinucleotide, Reduced Form, Preweighed Vial, Disodium Salt, Sigma Stock No. 340-110, in the appropriate volume of deionized water. **PREPARE FRESH.**)
- D. 1.0% (w/v) Bovine Serum Albumin (BSA)
(Prepare 10 ml in deionized water using Albumin, Bovine, Sigma Prod. No. A-6003.)
- E. Sorbitol Dehydrogenase Enzyme Solution
(Prepare a solution containing 70 - 150 units/ml of Sorbitol Dehydrogenase in cold deionized water. Store at 4°C for 1 hour. Immediately before use, dilute to a final concentration of 0.55 - 0.75 unit/ml with cold Reagent D.)

**Enzymatic Assay of SORBITOL DEHYDROGENASE
(EC 1.1.1.14)**

PROCEDURE:

Pipette (in milliliters) the following reagents into suitable cuvettes:

	<u>Test</u>	<u>Blank</u>
Reagent A (Buffer)	2.35	2.35
Reagent B (Fructose)	0.50	0.50
Reagent C (β -NADH)	0.05	0.05

Mix by inversion and equilibrate to 25°C. Monitor the A_{340nm} until constant, using a suitably thermostatted spectrophotometer. Then add:

Reagent E (Enzyme Solution)	0.10	-----
Reagent D (BSA)	-----	0.10

Immediately mix by inversion and record the decrease in A_{340nm} for approximately 5 minutes. Obtain the $\Delta A_{340nm}/\text{minute}$ using the maximum linear rate for both the Test and Blank.

CALCULATIONS:

$$\text{Units/ml enzyme} = \frac{(\Delta A_{340nm}/\text{min Test} - \Delta A_{340nm}/\text{min Blank})(3)(df)}{(6.22)(0.1)}$$

3 = Total volume (in milliliters) of assay

df = Dilution factor

6.22 = Millimolar extinction coefficient of β -NADH at 340nm

0.1 = Volume (in milliliter) of enzyme

$$\text{Units/mg solid} = \frac{\text{units/ml enzyme}}{\text{mg solid/ml enzyme}}$$

$$\text{Units/mg protein} = \frac{\text{units/ml enzyme}}{\text{mg protein/ml enzyme}}$$

UNIT DEFINITION:

One unit will convert 1.0 μmole of D-fructose to D-sorbitol per minute at pH 7.6 at 25°C.

**Enzymatic Assay of SORBITOL DEHYDROGENASE
(EC 1.1.1.14)**

FINAL ASSAY CONCENTRATION:

In a 3.00 ml reaction mix, the final concentrations are 78 mM triethanolamine, 183 mM D-fructose, 0.2 mM β -nicotinamide adenine dinucleotide, reduced form, 0.033% (w/v) bovine serum albumin and 0.055 - 0.075 unit sorbitol dehydrogenase.

REFERENCE:

Gerlach, U. and Hiby, W. (1974) in *Methods of Enzymatic Analysis* (Bergmeyer, H.U. ed.) 2nd ed., Volume II, 569-573, Academic Press Inc., New York, NY

NOTES:

1. This assay is based on the cited reference.
2. Where Sigma Product or Stock numbers are specified, equivalent reagents may be substituted.

Sigma warrants that the above procedure information is currently utilized at Sigma and that all Sigma-Aldrich, Inc. products conform to the information in this and other Sigma-Aldrich, Inc. publications. Purchaser must determine the suitability of the information and product(s) for their particular use. Additional terms and conditions may apply. Please see reverse side of the invoice or packing slip.