

Quick Start Guide

mPAGE® TurboMix Bis-Tris Gel Casting Kit

TMKIT-10
TMKIT-60



5 Easy Steps to Cast

The mPAGE® TurboMix Bis-Tris Gel Casting Kit uses a quick cast method which can be used to prepare protein separation gels for SDS-PAGE. The volumes provided in this protocol are for casting a single 7.4 x 8.2 cm mini gel, for different gel sizes solution volumes must be determined by the user.


For additional gel sizes and traditional casting instructions, read the complete User Guide on the mPAGE® TurboMix Bis-tris Gel Casting Kit product page at [SigmaAldrich.com](https://www.sigmaaldrich.com).

Read the complete user guide before first use.

Bring reagents to room temperature prior to use.
 Important: Bis-Tris gels are only compatible with MOPS-SDS and MES-SDS running buffer. Do NOT use Tris-Glycine running buffer.

1. Prepare Resolving Gel


Use the table below to determine solution volumes. Gently mix all reagents in a clean 15 mL conical tube.

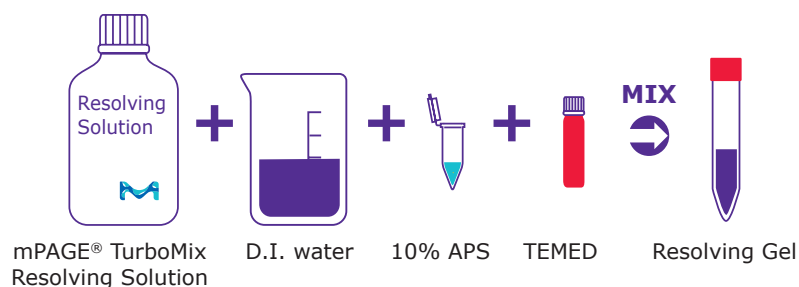

Note: Gel will begin to polymerize after addition of APS and TEMED.

2. Prepare Stacking Gel

Gently mix all reagents in a clean 15 mL conical tube.

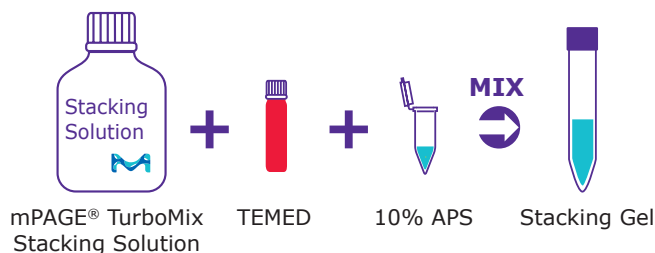

Volumes required to cast one 7.4 x 8.2 cm mini gel using the mPAGE® Gel Caster:

- For 0.75 mm gels, multiply volumes in table by 0.75.
- For 1.50 mm gels, multiply volumes in table by 1.5.

Resolving Gel

Mini Gel Thickness 1 mm


Gel percentage	8%	10%	12%	15%
mPage® TurboMix Resolving Solution	2.4 mL	3 mL	3.6 mL	4.5 mL
D.I. water	3.6 mL	3 mL	2.4 mL	1.5 mL
10% APS	30 µL	30 µL	30 µL	30 µL
TEMED	3 µL	3 µL	3 µL	3 µL
Total	6 mL	6 mL	6 mL	6 mL

Stacking Gel

Mini Gel Thickness 1 mm


Gel percentage	4%
mPage® TurboMix Stacking Solution	2 mL
D.I. water	--
10% APS	20 µL
TEMED	2 µL
Total	2 mL

3. Add Resolving Gel

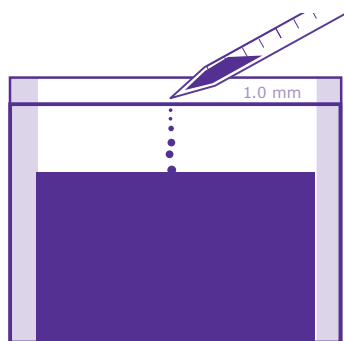


Using a serological pipette, fill casting cassette with resolving gel to desired height.

4. Add Stacking Gel

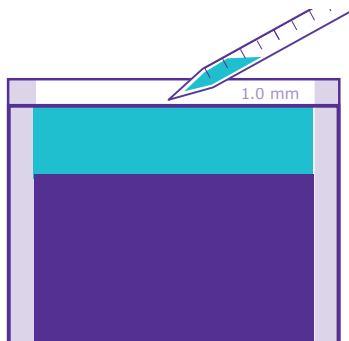

Using a serological pipette, slowly add stacking gel solution until reaching top of the shorter glass plate.

5. Allow to Polymerize

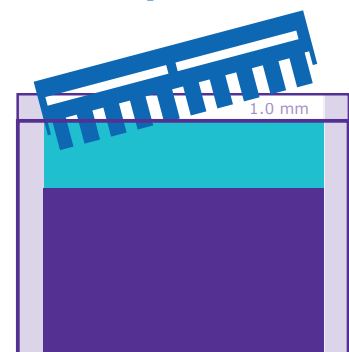

Insert comb and allow gel to polymerize, 1 hour.

Notice

We provide information and advice to our customers on application technologies and regulatory matters to the best of our knowledge and ability, but without obligation or liability. Existing laws and regulations are to be observed in all cases by our customers. This also applies in respect to any rights of third parties. Our information and advice do not relieve our customers of their own responsibility for checking the suitability of our products for the envisaged purpose.

The information in this document is subject to change without notice and should not be construed as a commitment by the manufacturing or selling entity, or an affiliate. We assume no responsibility for any errors that may appear in this document.

Technical Assistance

Visit the tech service page on our web site at SigmaAldrich.com/TechService.

Terms and Conditions of Sale

Warranty, use restrictions, and other conditions of sale may be found at SigmaAldrich.com/Terms.

Contact Information

For the location of the office nearest you, go to SigmaAldrich.com/Offices.

MilliporeSigma, Millipore, mPAGE and Sigma-Aldrich are trademarks of Merck KGaA, Darmstadt, Germany or its affiliates. All other trademarks are the property of their respective owners. Detailed information on trademarks is available via publicly accessible resources.
© 2022 Merck KGaA, Darmstadt, Germany and/or its affiliates. All Rights Reserved.

The life science business of Merck KGaA, Darmstadt, Germany operates as MilliporeSigma in the U.S. and Canada.