

## Technical Data Sheet

# Tetrathionate-Brilliant-Green (TBG) Bile Enrichment Broth, modified for microbiology

Ordering number: 1.05178.0500

Tetrathionate-brilliant-green (TBG) bile enrichment broth is used for the selective enrichment of *Salmonella* for the examination of pharmaceutical products in raw materials as well as foodstuffs, meat etc.

The medium complies with the recommendations of the APHA for the examination of food (1992).

### Mode of Action

Brilliant green, ox bile and high concentrations of thiosulfate and citrate largely inhibit the accompanying microbial flora. Sulfide production is detected by using thiosulfate and iron ions, the colonies turn black. The presence of *coliform* bacteria is established by detecting degradation of lactose to acid with the pH indicator neutral red.

### Typical Composition

Peptones	10 g/l
Lactose	10 g/l
Ox Bile	8.5 g/l
Sodium Citrate	10 g/l
Na <sub>2</sub> S <sub>2</sub> O <sub>3</sub>	8.5 g/l
Ammonium Iron(III) Citrate	1.0 g/l
Brilliant Green	0.0003 g/l
Neutral Red	0.025 g/l
Agar-Agar	12 g/l

### Preparation

Suspend 60 g/l completely. Pour plates. **Do not autoclave.**

The appearance of the plates is turbid with sediment and green.

The pH value at 25 °C is in the range of 6.8-7.2.

## Specimen

e.g. Stool,

Clinical specimen collection, handling and processing, see general instructions of use.

## Experimental Procedure and Evaluation

Spread the sample or material from an enrichment culture on the surface of the culture medium.

Incubation 24 h at 35 °C aerobically.

Lactose-negative colonies are colorless. Lactose-positive colonies are pink to red. Colonies of microorganisms producing H<sub>2</sub>S have a black center.

Appearance of Colonies	Microorganisms
Colourless, translucent	<i>Shigella</i> and some <i>Salmonella</i> species
Translucent with a black centre	<i>Proteus</i> and most <i>Salmonella</i> species
Pink to red	<i>Escherichia coli</i>
Colonies are larger than those of <i>Escherichia coli</i> , pink to whitish or cream-coloured, opaque, mucoid	<i>Enterobacter aerogenes</i>

## Storage

The product can be used for sampling until the expiry date if stored upright, protected from light and properly sealed at +15 °C to +25 °C.

After first opening of the bottle the content can be used up to the expiry date when stored dry and tightly closed at +15 to +25° C.

## Disposal

Please mind the respective regulations for the disposal of used culture medium (e.g. autoclave for 20 min at 121 °C, disinfect, incinerate etc.).

## Quality Control

Control Strains	ATCC #	Inoculum CFU	Incubation	Expected Results
<i>Salmonella typhimurium</i>	14028	10-100	24 h at 35 °C	Enrichment from mixed inoculum ≥ 95 %, growth
<i>Escherichia coli</i>	25922	≥ 10 <sup>4</sup>	24 h at 35 °C	Enrichment from mixed inoculum ≤ 5 %, growth inhibited

Please refer to the actual batch related Certificate of Analysis.



We provide information and advice to our customers on application technologies and regulatory matters to the best of our knowledge and ability, but without obligation or liability. Existing laws and regulations are to be observed in all cases by our customers. This also applies in respect to any rights of third parties. Our information and advice do not relieve our customers of their own responsibility for checking the suitability of our products for the envisaged purpose.

EMD Millipore and the M logo are registered trademarks of Merck KGaA, Darmstadt, Germany. ATCC is a registered trademark of ATCC, Manassas, VA, USA. Lit. No. TN7413ENUS

## Literature

American Public Health Association (1992). Compendium of methods for the microbiological examination of foods. 3<sup>rd</sup> edition.

Deutsches Arzneibuch (DAB), 10. Auflage, Kapitel VIII, 10.

European Pharmacopeia II (2003). Chapter VIII, 10.

Jeffries, L. (1959). Novobiocin-tetrathionate broth: A medium of improved selectivity for the isolation of *salmonellae* from faeces. J. Clin. Path., **12**: 568-571.

## Ordering Information

Product	Cat. No.	Pack size
TBG-Broth modified for microbiology (Tetra-thionate-Brilliant-green Bile Enrichment Broth)	1.05178.0500	500 g
BPLS Agar for the isolation of <i>Salmonella</i>	1.07237.0500	500 g
Lactose broth for microbiology	1.07661.0500	500 g
LEIFSON Agar (Deoxycholate citrate agar acc. to LEIFSON, modified for microbiology)	1.02896.0500	500 g
XLD Agar (acc. harm. EP/USP/JP)	1.05290.0500	500 g

EMD Millipore Corporation  
290 Concord Road  
Billerica, MA 01821, U.S.A.  
mibio@emdmillipore.com  
[www.emdmillipore.com/biomonitoring](http://www.emdmillipore.com/biomonitoring)

Find contact information for your country at:  
[www.emdmillipore.com/offices](http://www.emdmillipore.com/offices)  
For Technical Service, please visit:  
[www.emdmillipore.com/techservice](http://www.emdmillipore.com/techservice)



We provide information and advice to our customers on application technologies and regulatory matters to the best of our knowledge and ability, but without obligation or liability. Existing laws and regulations are to be observed in all cases by our customers. This also applies in respect to any rights of third parties. Our information and advice do not relieve our customers of their own responsibility for checking the suitability of our products for the envisaged purpose.

EMD Millipore and the M logo are registered trademarks of Merck KGaA, Darmstadt, Germany. ATCC is a registered trademark of ATCC, Manassas, VA, USA. Lit. No. TN7413ENUS