

Product Information

Carbonic Anhydrase II, human recombinant, expressed in *Escherichia coli*

Catalog Number **C6624**
Storage Temperature $-20\text{ }^{\circ}\text{C}$

Synonym: CA II

Product Description

Carbonic anhydrases (CA) are a family of enzymes that catalyze the rapid conversion of carbon dioxide to bicarbonate and protons, a reaction that occurs rather slowly in the absence of a catalyst. The active site of most carbonic anhydrases contains a zinc ion. They are, therefore, classified as metalloenzymes.¹⁻²

Carbonic anhydrases are widely distributed in plant and animal tissues where they are involved in diverse physiological processes, such as photosynthesis, pH homeostasis, calcification, and bone resorption.

There are at least five distinct CA families (α , β , γ , δ , and ϵ). These families have no significant amino acid sequence similarity and, in most cases, are thought to be an example of convergent evolution. The α -CAs are found in humans. At least 14 isoforms of α -CA have been identified in vertebrates with different physiological and pathological roles. The enzymes can be localized in the cytosol or mitochondria, be membrane bound with extracellular domains, or be secreted.²⁻⁴


Carbonic anhydrase II (CA II) is a cytosolic, ~29 kDa protein. It is required for kidney renal acidification and its absence in an inherited syndrome leads to osteoporosis, renal tubular acidosis, and cerebral calcification.⁵⁻⁷ In addition, it has long been a drug target for the treatment of glaucoma, but has taken on new importance as potential target for obesity and cancer.⁸⁻¹²

The product is supplied as a solution in 20 mM Tris, pH 7.5, with 150 mM NaCl.

Purity: $\geq 90\%$ (SDS-PAGE)

Enzyme activity:


1. Conversion of carbon dioxide to bicarbonate:

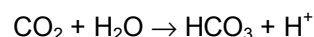

Specific activity: $\geq 5,000$ units/mg

Unit Definition: One unit will decrease the pH of a 20 mM Tris buffer from pH 8.3 to 6.3 in 1 minute at $0\text{ }^{\circ}\text{C}$.

2. Esterase activity:

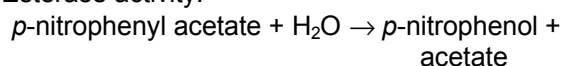

The increase in the amount of the product, *p*-nitrophenol, is measured as absorbance increases at 405 nm.

Specific activity: ≥ 2 nmole/min/ μg

Precautions and Disclaimer

This product is for R&D use only, not for drug, household, or other uses. Please consult the Material Safety Data Sheet for information regarding hazards and safe handling practices.

Storage/Stability

Store the product at $-20\text{ }^{\circ}\text{C}$. The product is stable for at least 2 years as supplied.

After initial thawing, the enzyme should be refrozen at $-20\text{ }^{\circ}\text{C}$ in aliquots.

References

1. Badger, M.R., and Price, G.D., The role of carbonic anhydrase in photosynthesis. *Annu. Rev. Plant Physiol. Plant Mol. Bio.*, **45**, 369-392 (1994).
2. Tashian, R.E., The carbonic anhydrases: widening perspectives on their evolution, expression and function. *BioEssays*, **10**, 186-192 (1989).
3. Tashian, R.E., Genetics of the mammalian carbonic anhydrases. *Adv. Genet.*, **30**, 321-356 (1992).

4. Kavroulakis, N., et al., Tissue distribution and subcellular localization of carbonic anhydrase in mature soybean root nodules indicates a role in CO₂ diffusion. *Plant Physiol. Biochem.*, **41**, 479-484 (2003).
5. Avvaru, B.S., et al., Apo-Human carbonic anhydrase II revisited: implications of the loss of a metal in protein structure, stability, and solvent network. *Biochemistry*, **48**, 7365-7372 (2009).
6. Borthwick, K.J., et al., A phenocopy of CAII deficiency: a novel genetic explanation for inherited infantile osteopetrosis with distal renal tubular acidosis. *J. Med. Genet.*, **40**, 115-121 (2003).
7. Whyte, M.P., Carbonic Anhydrase II Deficiency. *Clin. Orthop. Relat. Res.*, **294**, 52-63 (1993).
8. Recacha, R., et al., Crystal structure of human carbonic anhydrase II complexed with an anti-convulsant sugar sulphamate. *Biochem. J.*, **361**, 437-441 (2002).
9. Casini, A., et al., Carbonic anhydrase inhibitors: topically acting antiglaucoma sulfonamides incorporating esters and amides of 3- and 4-carboxybenzamide. *Bioorg. Med. Chem. Lett.*, **13**, 2867-2873 (2003).
10. Supuran, C.T., Carbonic anhydrase inhibitors in the treatment and prophylaxis of obesity. *Expert. Opin. Ther. Pa.*, **13**, 1545-1550 (2003).
11. Ho, Y.T., et al., Inhibition of carbonic anhydrase II by steroidal and non-steroidal sulphamates. *Biochem. Biophys. Res. Commun.*, **305**, 909-914 (2003).
12. Abbate, F., et al., Carbonic anhydrase inhibitors: X-ray crystallographic structure of the adduct of human isozyme II with EMATE, a dual inhibitor of carbonic anhydrase and steroid sulfatase. *Bioorg. Med. Chem. Lett.*, **14**, 231-234 (2004).

EB, RBG, MAM 08/10-1

Sigma brand products are sold through Sigma-Aldrich, Inc.

Sigma-Aldrich, Inc. warrants that its products conform to the information contained in this and other Sigma-Aldrich publications. Purchaser must determine the suitability of the product(s) for their particular use. Additional terms and conditions may apply. Please see reverse side of the invoice or packing slip.